

Separation of Selected Components of the Allergens Contained in Balsam of Peru

The Essential Oil Balsam of Peru is commonly found in a multitude of products, including, but not limited to: perfumes, cosmetics, sunscreens, and flavored beverages. The fragrant liquid is harvested from trees grown in Central America and smells richly of cinnamon, vanilla, and clove. The oil has been used in three main areas: as a medicinal product pertaining to its anti-septic, anti-fungal and anti-parasitic natures, as a flavor enhancer for food and drinks, and as a fragrance in perfumes and toiletries. Despite the utility of the mixture, it has been shown that allergies can arise from some of the individual components. Common symptoms generally present in the form of contact dermatitis and eczema. Use of such components in the manufacturing of the above mentioned products requires quantification of the individual compounds. Therefore, we have developed an HPLC method that separates five components known to have allergenic properties. The method can be used with either the PRP-C18 (5 µm) or the PRP-1 (5 µm) column. The two chromatograms illustrate the versatility and selectivity of the PS-DVB backbone found in all of Hamilton Company columns.

Column Information


Packing Material	Dimensions	Part Number
PRP-1 (5 µm)	150 x 4.6 mm	79641
PRP-C18 (5 µm)	150 x 4.6 mm	79676

Chromatographic Conditions

Gradient	0.00–2 min, 75% B
	2.01–5 min, 75–100% B
	5.01–8 min, 100% B
Temperature	Ambient
Injection Volume	10 µL
Detection	UV at 214 nm
Eluent A	Water
Eluent B	Acetonitrile
Flow Rate	2.0 mL/min (PRP-1)
	2.5 mL/min (PRP-C18)

Author: Adam L. Moore, PhD, Hamilton Company


PRP-1 Separation of Esters in Balsam of Peru

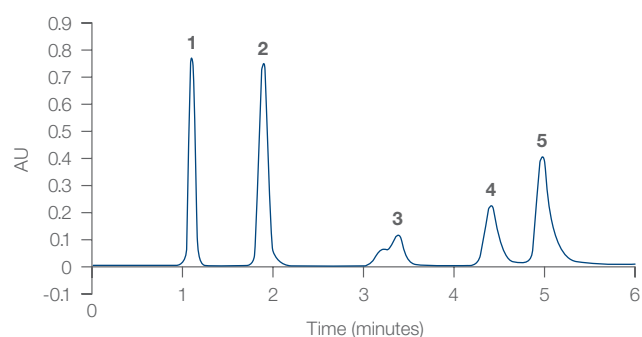

Compounds:

1: Benzyl Alcohol 3: Farnesol 5: Benzyl Cinnamate
2: Trans-Cinnamaldehyde 4: Benzyl Benzoate

PRP-C18 Separation of Esters in Balsam of Peru

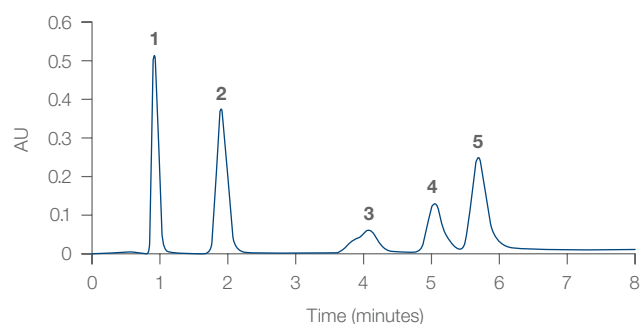

Compounds:

1: Benzyl Alcohol 3: Farnesol 5: Benzyl Cinnamate
2: Trans-Cinnamaldehyde 4: Benzyl Benzoate

©2019 Hamilton Company. All rights reserved.
All other trademarks are owned and/or registered by Hamilton Company in the U.S. and/or other countries.
Lit. No. L80100 – 09/2019

Web: www.hamiltoncompany.com

USA: 800-648-5950

Europe: +40-356-635-055

To find a representative in your area, please visit hamiltoncompany.com/contacts.

Hamilton Americas & Pacific Rim

Hamilton Company Inc.
4970 Energy Way
Reno, Nevada 89502 USA
Tel: +1-775-858-3000
Fax: +1-775-856-7259
sales@hamiltoncompany.com

Hamilton Europe, Asia & Africa

Hamilton Central Europe S.R.L.
str. Hamilton no. 2-4
307210 Giarmata, Romania
Tel: +40-356-635-055
Fax: +40-356-635-060
contact.lab.ro@hamilton-ce.com