

Introduction

Key benefits of SGE syringes are:

- Longer life
- Superior performance and robustness
- Reduced carryover

In order to maintain optimum performance and long life, a few simple care issues should be considered.

Instructions

- For optimal performance, it is recommended that eVol® syringes only be used in the temperature range of 21 - 23°C.
- Always inspect the syringe before use. Check the barrel for cracks and the needle tip for burrs.
- To eliminate carryover between samples, flush the syringe with solvent 5-20 times, remembering to discard the first 2-3 washes.

Installing the syringe

- Pull up on the bottom collar of the eVol and insert the back (plunger end) of the syringe into the bottom of the instrument. Ensure the syringe plunger is in the zero position.
- If the syringe has not been used with the eVol unit, it will automatically re-zero the syringe.


Removing the syringe

- Pull up on the bottom collar of the eVol to remove the syringe.
- Pull the syringe out. The plunger will be extended as you pull out the syringe and it disengages from the magnet.


Clean and care

- Syringes should be cleaned thoroughly after use.
- Syringe cleaning agents will usually depend on the contaminating material. Methanol, methylene chloride, acetonitrile and acetone are commonly used solvents for cleaning.
- Clean externally by wiping with a tissue.

- After the syringe has been cleaned, rinse the syringe with acetone, remove plunger and air dry.

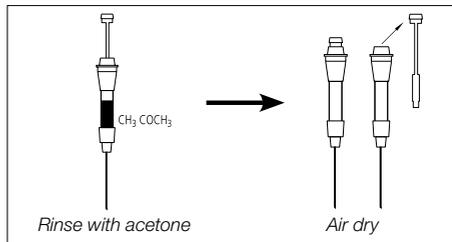

- Sterilization modes:
 - Autoclave: Remove plunger from the barrel prior to autoclaving. Set autoclave to a maximum temperature of 100°C. Allow syringes to come to room temperature before inserting the plunger back into the barrel.
 - Gas: Use ethylene oxide
- Do not immerse the entire syringe in solvent as this may damage the adhesive used to bond parts of the syringe.

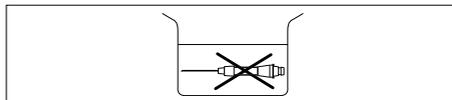


Plunger care

- Avoid cycling the plunger when the syringe is dry.
- To remove plunger unscrew black plunger guide.

Needle care

- Check the needle for burrs prior to use. Use a fine emery board or carborundum to remove burrs.
- To unblock needles, remove the plunger and fill the syringe with solvent using another syringe. Re-insert the plunger and gently push solvent through needle. Never force the plunger. Too much pressure may crack the syringe barrel.

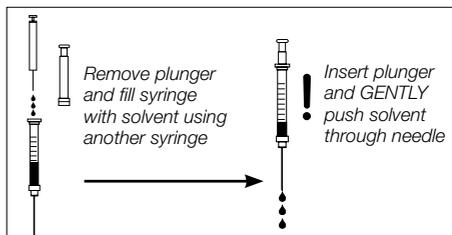

- The following syringe needle cleaning product is available:
 - Needle cleaning kit: The kit contains a range of stylet wires for needle cleaning, tweezers and a surfactant material for barrel cleaning.

Replacement

Replacement needles

- To replace needles unscrew front cover nut and spring, remove needle. Carefully insert replacement needle into the front of the syringe, slide spring and cover nut over the needle and screw nut onto the front screw of the syringe barrel.
- For replacement needles refer to www.trajanscimed.com

Storage

- To protect against breakage store the syringe in its original packaging or on a syringe rack.
- Always thoroughly clean the syringe and ensure that it is dry before storage.

Warranty

All Trajan Scientific and Medical syringes are warranted to meet the stated quality and performance specifications and to be free of defects in material or workmanship. The warranty implies free replacement of a defective syringe only upon proper written proof of the defect and if requested by Trajan Scientific and Medical the return of the defective product in its original packaging. It does not apply to mishandling of product by the customer, either in storage or use.

- Do not return any syringe that has been used with radioactive, infectious or hazardous materials.
- Do not return any syringe unless requested to do so by an authorized Trajan representative.
- Any syringe being returned must be clearly marked with a return authorization number, which is available from your Trajan representative.

Conformance statement

Trajan syringes are manufactured under a documented Quality Management System. All in-process measuring equipment and instrumentation is maintained and calibrated in accordance with stringent quality standards. Trajan warrants syringe displacement within the stated conformity specifications.

Information and support

Visit www.trajanscimed.com or contact techsupport@trajanscimed.com

Specifications are subject to change without notice.